

- 11. Update the Software
  - a. Plug in USB flash drive to the USB port located on the top, near the power inlet.



b. Power on the unit, then go to **Menu**.









c. Log in with **service** user account

Sa ( Pr	service	
	Your full name	
	Password	

d. Go to Service menu, then Software Update



e. Tap **Next** button. It will show the current firmware name and version. The below is only illustration. The actual current firmware name and version may be different.









f. Tap **OK** 



g. Wait for the updater page to open



h. Select Update via USB









i. Select the provided update file. The update file shall have extention \*.swu .

Update via USB		
Selected software:	file:///D:/AhmaupdateApp/debug	
	Up Support: FAT16, FAT32 and ExFAT	
General Back	Update	

j. Then the selected update file will be displayed on the left side. Make sure the software name is the same with the existing software name but different version number. So that you will not require to do any recertification process after a software update.

	Update vi	a USB
	Selected software: Esco-HMI-Software-Update.swu	file:///D:/AhmaupdateApp/debug   Esco-HMI-Software-Update.swu  Up  Up  Support: FAT16, FAT32 and ExFAT
€	Back	Update

1. Tap Update

k.









m. Wait for the update process to finish



n. When the update is finish, unplug the USB flash drive and press Restart.





